

透明化试剂选择流程

1. 按目标组织选择

包括脑在内的所有组织：CUBIC
 脑组织：SCALEVIEW-S, SeeDB 或 SeeDB2
 多细胞球状体：SCALEVIEW-S4
 植物：ClearSee

2. 按样品大小选择

全脑：CUBIC
 1mm 脑切片~半脑：SCALEVIEW-S
 0.5mm~2mm 脑切片：SeeDB 或 SeeDB2

3. 按分析方法选择

综合分析（激光片层扫描显微镜）：CUBIC
 高分辨率分析（双光子共聚焦显微镜）：SCALEVIEW-S, SeeDB 或 SeeDB2
 超高分辨率分析（超分辨显微镜）：SeeDB2

4. 按样品观察时间选择

长期（大于三周）：SCALEVIEW-S 或 SeeDB2
 中期（一至三周）：CUBIC（荧光蛋白信号会减弱）
 短期（一周）：SeeDB（在处理的一周后会出现自体荧光）

5. 按抗体染色选择

SCALEVIEW-S > CUBIC = SeeDB2

6. 按透明度选择

CUBIC > SCALEVIEW-S > SeeDB2 或 SeeDB

	CUBIC	SCALEVIEW-S	SeeDB	SeeDB2	ClearSee
组织	所有组织	脑组织, 球体组织	脑组织	脑组织	植物组织
样品大小	全脑	脑切片(1mm)~半脑	脑切片(0.5~2.0mm)		叶, 茎, 根等
分析方法	综合分析 (激光片层扫描显微镜)	高分辨率分析(双光子共聚焦显微镜)			
时间长短	中期(1-3周)	长期(3周以上)	短期(1周)	长期(3周以上)	长期(3周以上)

上述试剂仅供实验研究用,不可用作“医药品”、“食品”、“临床诊断”等。

Listed products are intended for laboratory research use only, and not to be used for drug, food or human use. / Please visit our online catalog to search for other products from FUJIFILM Wako; <https://labchem-wako.fujifilm.com/> / This leaflet may contain products that cannot be exported to your country due to regulations. / Bulk quote requests for some products are welcomed. Please contact us.

富士胶片和光(广州)贸易有限公司 labchem.fujifilm-wako.com.cn
 广州市越秀区先烈中路69号东山广场30楼 3002-3003室
 北京 Tel: 13611333218
 上海 Tel: 021 62884751
 广州 Tel: 020 87326381
 香港 Tel: 852 27999019
 询价: wkgz.info@fujifilm.com
 目录价查询: bb-china.net

官方微信



目录价查询

